

ES

Provided with an Automatic Air Release Time Setting Function

HIGH PRESSURE STEAM STERILIZER ES-315 / ES-215

Reliable & Speedy Sterilization

Automatic air release time setting function allows an optimal evacuation time for the remaining air in the chamber to be set automatically according to the type and quantity of article to be sterilized, and ensures more reliable and speedy sterilization. This function is equipped with ES-Series as a standard feature.

User-friendly Memory Function

Once turned on the system automatically sets and displays a pre-set value. Operation is done simply by pressing the start switch. ES-315/ES-215 enables the set temperature and time for each mode program to be stored in its memory.

Easy-to-see Control Panel



※The control panel lighting image shown above may vary from actual display.

MODE/PROCESS display enables real time monitoring of operation status.
————— ES-315 / ES-215 offers three mode programs. —————


Sterilization Mode

It contains a delay start timer function that allows operation to start 1 to 99 hours later.



Heating Mode

This mode is useful when not removing culture media from the autoclave immediately after sterilization. Sterilized media can be kept warm to prevent coagulation.

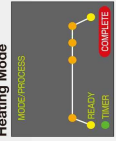

This mode program automates the dissolution of culture media that can sometimes be a painful process when using a water bath to boil them.

User-friendly design allows for even more easy maintenance.

Chamber having a flat inner surface

All models are supplied with stainless steel wire mesh baskets. No protrusions on inner surface of the chamber enable smooth loading and unloading of the basket as well as block design of all pipe fittings ensures safe operation.

Front access for easy maintenance

The exhaust bottle to capture and store exhaust steam is placed on the front panel of the autoclave unit. Access to the chamber for replacing water inside, as well as for routine maintenance and inspection, is done on the front of the unit, which also permits operators to replace bellows, saving costs for services by the service personnel.

A microprocessor controlled monitoring system provides safe sterilization.

Incomplete sterilization protection

In case of abnormal chamber temperature rise, the autoclave stops its operation. In case of abnormal decrease in temperature, the sterilization timer is temporarily suspended and then reactivated when the set temperature is reached.

Chamber overpressure protection

If overpressure in the chamber is detected, the autoclave stops all the function and indicates an error message simultaneously.

Chamber interior overtemperature protection

If the temperature inside the chamber exceeds the set temperature by 5 °C or more, the autoclave stops all the function and indicates an error message simultaneously to avoid overheating.

Chamber exterior overtemperature protection


The thermostat switch detects the temperature of chamber exterior wall, which allows the autoclave to stop all the function if the temperature exceeds 130°C.

◆ Specifications

Model	ES-215	ES-315
Operating temperature range (during sterilizing)	105 - 123 °C	105 - 132 °C
Operating pressure range	0 - 127kPa	0 - 188kPa
Maximum operating pressure	147kPa	216kPa
Temperature control	Digital, microprocessor controlled	Digital, microprocessor controlled
Temperature display / display range	Digital / -15 - 180 °C	Digital / -15 - 180 °C
Pressure display / display range	Analog / 0 - 250kPa	Analog / 0 - 400kPa
Heat source	1.5kW electric heater	2.0kW electric heater
Timer control	Digital, microprocessor controlled	Digital, microprocessor controlled
Timer display	Digital	Digital
Timer display range	1 - 240min. (sterilizing) / 1 - 8hours (warming)	1 - 240min. (sterilizing) / 1 - 8hours (warming)
Start-timer setting range	1 - 99hours in each operation mode	1 - 99hours in each operation mode
Memory function	A set of temperature and time setting in each operation mode	A set of temperature and time setting in each operation mode
Safety devices	● Inside the chamber overheat prevention, ● Outer wall of the chamber overheat prevention, ● Overpressure prevention, ● Temperature sensor disconnection prevention, ● Empty heating prevention, ● Leakage breaker, ● Safety valve	● Inside the chamber overheat prevention, ● Outer wall of the chamber overheat prevention, ● Overpressure prevention, ● Temperature sensor disconnection prevention, ● Empty heating prevention, ● Leakage breaker, ● Safety valve
Malfunction prevention devices	● Low water level detection, ● Exhaust valve knob open/close detection, ● Insufficient sterilization detection, ● Chamber lid open/close detection	● Low water level detection, ● Exhaust valve knob open/close detection, ● Insufficient sterilization detection, ● Chamber lid open/close detection
Leakage breaker	20A (120V) , 15A (220 / 230 / 240V)	30A (120V) , 15A (220 / 230 / 240V)
Rated breaking current:	30mA (120V) , 10mA (220 / 230 / 240V)	30mA (120V) , 10mA (220 / 230 / 240V)
Rated sensed current:	Class 1 equipment	Class 1 equipment
Protection type	5 - 35 °C	5 - 35 °C
Operating environment	Relative humidity: 30 - 85%	Relative humidity: 30 - 85%
Atmospheric pressure:	700 - 1060hPa	700 - 1060hPa
Gradient:	Within 2°	Within 2°
Dimensions	φ248 x 543mm	φ325 x 740mm
Usable dimension	φ238 x 450mm	φ312 x 635mm
Capacity of chamber	22 ℓ	53 ℓ
Chamber material	SUS304	SUS304
Capacity / material of exhaust bottle	3 ℓ / polyethylene	3 ℓ / polyethylene
Type / material of lid gasket	Accommodating internal pressure / silicone rubber	Accommodating internal pressure / silicone rubber
Dimensions of main unit (mm)	400W x 460D x 920H (Height from floor to control panel/705)	490W x 960D x 1050H (Height from floor to control panel/875)
Net weight	50kg	80kg
Rated Voltage	AC 120 / 220 / 240V	AC 120 / 220 / 240V
Power input	15A (120V), 7A (220 / 230 / 240V)	20A (120V), 10A (220V), 9A (230 / 240V)
Power supply requirements	Single-phase 120V AC (50 / 60Hz) 15A or above	Single-phase 120V AC (50 / 60Hz) 20A or above
Power Consumption (calorific value)	Single-phase 220 / 230 / 240V AC (50 / 60Hz) 15A or above	Single-phase 220 / 230 / 240V AC (50 / 60Hz) 15A or above
Accessories	1.5kW (120kcal/h) ● Stainless steel basket with slatted bottom plate 1, ● Water level sensor 1, ● Inspection sheet 1, ● Warranty card 1, ● Customer card 1, ● Operator's manual 1, ● Clear plastic accessory case 1, ● Attachment screw for clear plastic accessory case 1	2kW (1720kcal/h) ● Stainless steel basket with slatted bottom plate 1, ● Water level sensor 1, ● Inspection sheet 1, ● Warranty card 1, ● Customer card 1, ● Operator's manual 1, ● Clear plastic accessory case 1, ● Attachment screw for clear plastic accessory case 1

◆ For dimension details on the chamber and usable or inside dimension, please refer to each part dimensional drawing on the back page.

◆Part dimensional drawings

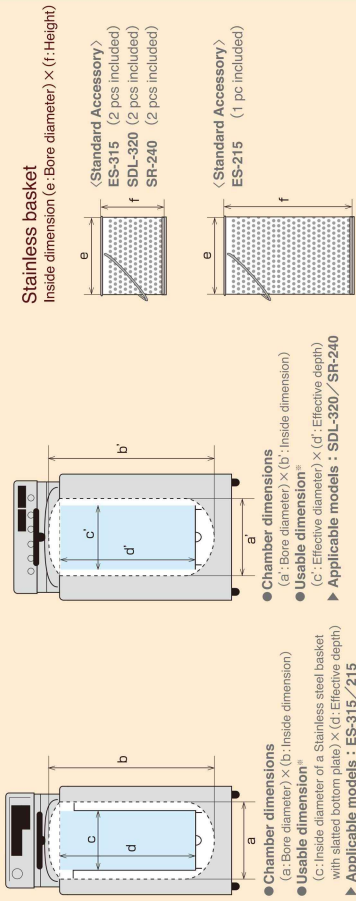

* The usable dimension shows recommended values considering safety during operation. The volume or position of the sterilization items may block the holes inside the chamber or interrupt the flow of steam or air, which will result in abnormal pressure increase or incomplete sterilization. To ensure safe operation of the autoclave, it is recommended that Stainless steel wire mesh baskets supplied with the unit as standard accessories or optional stainless steel buckets be used.

◆ The image shown for the product may not display its real color or shape.

◆Accessories



Stainless Basket

The entire basket is made of punched metal and is designed for easy penetration of steam.



Stainless Long Basket

Optimal for sterilizing the sterilizing bag. The structure of the lower section is without holes.



Stainless Bucket

Bucket without holes to prevent leakage of liquid from articles being sterilized.



Stand for Testing Durham's Tubes

Convenient for sterilizing test tubes containing culture media. Can be used as a stand independently.



Sterilizing Bag

Sterilizing bags fit to the vessel dimensions of each model are available.



Deodorant for autoclave

It allows for repeated use and removes offensive odors such as odors caused by culture media in the chamber.



Autoclavable waste container "ST-ZERO"

Waste inside the BSC can be collected in the container and autoclaved.

Sales Office

TOMY DIGITAL BIOLOGY CO., LTD.

3-14-17 Tagara, Nerima-ku, Tokyo 179-0073, Japan
e-mail : info@tomytech.com
URL : <http://www.digital-biology.co.jp>
phone : +81-3-5971-8160 fax : +81-3-3970-6036

TOMY TECH U.S.A., INC.

40479 Encyclopedia Circle, Fremont, California 94538, U.S.A.
e-mail : info@tomytech.com

URL : <http://www.tomytech.com>

phone : 510-440-1976 fax : 510-440-1975

Toll-Free US & Canada : 800-545-TOMT

TOMY SEIKO CO.,LTD.

Manufacturer

TOMY KOGYO CO., LTD.

3-14-17 Tagara, Nerima-ku, Tokyo 179-0073, Japan

All TOMY products have a limited one-year warranty. Specifications are subject to change according to product advancement. Tomy and Digital Biology is registered trademark of Tomy Seiko Co., Ltd. Tomy and Digital Biology are registered trademarks of Tomy Seiko Co., Ltd. Tomy Seiko and its subsidiaries. Printed in Japan.

3K08221N

TOMY



AUTOCLAVES

Sliding Door Type Series

HIGH PRESSURE STEAM STERILIZER
ES-315/215



SPEEDY AUTOCLAVE WITH DRIER
SDL-320



RETORT FOOD AUTOCLAVE
SR-240

